

# PCR Cabinet : UV Sterilisation

Sterilisation by UV radiation destroys unwanted PCR products

## Features

- Sterilisation by UV radiation
- Destroys unwanted PCR products
- Protects samples from contamination by user
- Built in timer control
- Doors are fitted with safety cutout switches
- Two powerful UV tubes provide efficient irradiation of the work surface

## Instructions for use

### Doors

Both doors must be closed and latched before UV light can be activated

### Timer for UV lights

Will run for up to 60 minutes. To set the timer, it must be turned fully to 60 then dialed back to the time required. If the timer is not fully turned before setting the time, the UV light will not run for the necessary time.

### Replacement tubes

In event of UV tube failure, please contact Thermoplastic Engineering for details of correct replacement tube.

## Specifications

Dimensions	W 600 x D 450 x H 755 mm
White light	8 watts
UV light	16 watts
Weight	15 kg
Voltage	230V AC

